

Series 4000 - Software and Data Removal

This document explains how to completely remove the Series 4000 software and data from your system. This might be necessary when encountering certain errors that prevent the software from starting up.

NOTE: All Environics data will be removed, including Cylinders, Concentration Mode settings, etc.

1. Close the Environics software
2. For Windows 7
 - a. Click **Start - Control Panel - Programs and Features**
 - b. Select Environics and click **Uninstall**
 - c. Open **Windows Explorer**. Click **Organize - Folder and Search Options**.
 - d. Click the **View** tab.
 - e. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.

For Windows 10

- a. Click **Start - Settings - Apps**
 - b. Select Environics and click **Uninstall**
 - c. Open **File Explorer**. Click **View - Options**.
 - d. Click the **View** tab.
 - e. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.
3. In Windows/File Explorer, navigate to the folder: **C:\programdata\environics**
4. Delete all files in folder: **C:\programdata\environics**

The Environics Software is now completely removed.