

## Series 4000 - Moving data file to another PC

This document explains how to move your Series 4000 data to a new PC, allowing you to keep all existing data

**IMPORTANT:** Be sure the same version of software is running on both the old and new PC, to avoid potential compatibility problems. If necessary, upgrade the software on the old PC first.

### ON OLD PC

1. Be sure the Environics software is closed.
2. For Windows 7
  - a. Open **Windows Explorer**. Click **Organize - Folder and Search Options**.
  - b. Click the **View** tab.
  - c. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.

#### For Windows 10

- a. Open **File Explorer**. Click **View - Options**.
  - b. Click the **View** tab.
  - c. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.
3. In Windows/File Explorer, navigate to the folder: **C:\programdata\environics**
4. Copy the file data.db to a USB thumb drive or Network drive

### ON NEW PC

1. Install the Environics software. Be sure the Environics software is not running.
2. For Windows 7
  - a. Open **Windows Explorer**. Click **Organize - Folder and Search Options**.
  - b. Click the **View** tab.
  - c. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.

#### For Windows 10

- a. Open **File Explorer**. Click **View - Options**.
  - b. Click the **View** tab.
  - c. Under **Advanced settings**, click **Show hidden files, folders, and drives**, and then click **OK**.
4. In Windows/File Explorer, navigate to the folder: **C:\programdata\environics**
5. Delete the file data.db
6. Copy the data.db file from the USB thumb drive to the C:\programdata\environics folder
7. Start the Environics software. All data from the old PC will be duplicated to the new PC